

Innovative Photonic Solutions, Inc. 4250 U.S. Highway 1, Suite 1 Monmouth Junction, NJ 08852 Phone: (732) 355-9300 Fax (732) 355-9302 sales@innovativephotonics.com http://www.innovativephotonics.com

Turn-key Integrated Raman Probe




Features

- 3-5X Higher Throughput than Standard Raman Probes (sample dependent)
- 785 nm Standard Wavelength Stabilized Excitation Source (514.5 nm, 830 nm, 1064 nm available)
- High Throughput Optical Design with 65cm⁻¹ Cut-on
- User Friendly Ergonomic Design
- Removable Distance Regulator¹ for Easy Sampling
- OEM Version Available

Innovative Photonic Solutions (IPS) is proud to introduce an ultra-high throughput Integrated Raman Probe. This novel device includes an integrated wavelength stabilized laser source with Raman filter packs, beam shaping optics and high efficiency Raman spectra collection optics. The probe interfaces with any fiber coupled spectrometer and simplifies operation and set-up.

The Integrated Raman Probe incorporates our wavelength stabilized hybrid external cavity laser (HECL) with a proprietary optical design to offer unmatched performance (typically 3 - 5x higher collection efficiency over traditional Raman probes). IPS's Integrated Raman Probe also comes complete with a UL/CE, and IEC compliant control box - providing a variety of power control options including modulation capability (TTL & analog) and a USB computer interface.

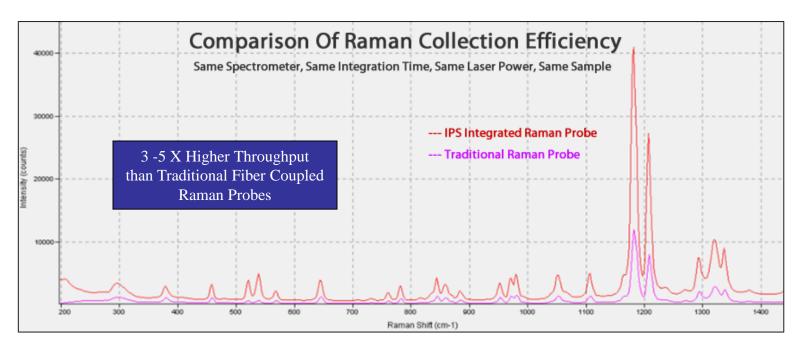

1 – Distance regulator available only for 9 mm working distance. 9 mm working distance is optimal if using IPS sample holder.

All data and statements contained herein are subject to change in accordance with Innovative Photonic Solution's policy of continual product improvement. No information contained herein is intended for use in connection with any contract except as may be first confirmed in writing by Innovative Photonic Solutions. The publication of information in this document does not imply freedom from patent or other rights of Innovative Photonic Solutions or others.



Innovative Photonic Solutions, Inc. 4250 U.S. Highway 1, Suite 1 Monmouth Junction, NJ 08852 Phone: (732) 355-9300 Fax (732) 355-9302 sales@innovativephotonics.com http://www.innovativephotonics.com


Integrated Raman Probe Technology & Specifications

Underlying Technology

IPS's Integrated Raman Probe offers higher collection efficiency as compared to traditional fiber probe approaches by optimizing the probe design in the following manner:

- IPS integrates the laser directly inside the probe head, eliminates fiber coupling losses and allows for beam shaping in order to optimize both laser power and power density on the sample which maximizes Raman signal.
- IPS utilizes a custom designed <u>rectangular core fiber</u> which increases both the coupling efficiency into both the fiber and the spectrometer.

Comparison of Loss at Fiber/Slit Interface & at Entrance to Collection Fiber (circle to line)

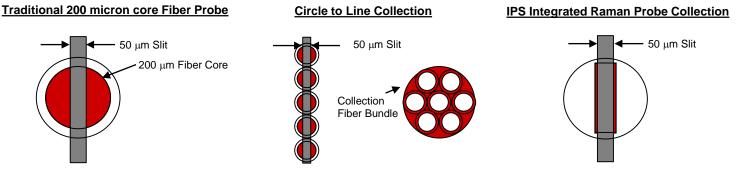

Raman signal is vignetted (thrown away) or not collected in red shaded regions detailed above

Parameter	Unit				
Excitation	Integral 785 nm wavelength stabilized laser <0.15 nm FWHM bandwidth				
	(514.5 nm, 830 nm and 1064 nm available)				
Collection	1.5 m long proprietary high throughput fiber				
Cut-on	65 cm-1 cut-on				
Electronic Connection	DB9 cable with safety interlock				
Power Control	Manual power adjustment knob, Analog / TTL modulation via BNC connector, or MicroUSB				
Power Supply	3 - 5 A max, 5VDC				
Shaft Material	316L Stainless Steel				
Fiber Bend Radius	4 inches				
Working distances	3 mm, 4 mm, 6.5 mm, and 9 mm standard (+/- 0.5 nm) - Custom distances available upon request				
Operating Temperature	15 degrees C to 35 degrees C				
Storage Temperature	- 20 degrees C to + 80 degrees C				
Humidity	0 - 80% non-condensing				

Parameter	Unit	Min	Тур	Max	Notes
Output power stability	%		± 1		
Peak wavelength	nm	784.5	785	785.5	
3 dB bandwidth (FWHM)	nm		0.1	0.15	
Operating Temperature Range (Case)	Deg C	15		35	Case Temperature
Power Consumption	W		3	7	Case temp between 15 and 35 deg C
Wavelength Stability	Seconds			180	Cold Start - to < 1 wavenumber
				1	Warm Start - to < 1 wavenumber
				3	Warm Start - to < 0.1 wavenumber

All data and statements contained herein are subject to change in accordance with Innovative Photonic Solution's policy of continual product improvement. No information contained herein is intended for use in connection with any contract except as may be first confirmed in writing by Innovative Photonic Solutions. The publication of information in this document does not imply freedom from patent or other rights of Innovative Photonic Solutions or others.




Innovative Photonic Solutions, Inc. 4250 U.S. Highway 1, Suite 1 Monmouth Junction, NJ 08852 Phone: (732) 355-9300 Fax (732) 355-9302 sales@innovativephotonics.com http://www.innovativephotonics.com

Standard 785 nm Integrated Raman Probes

Laser Type	Connector	Working Distance	Part Number
Single-Mode	SMA	3 mm	10785SP100-T030S
		4 mm	10785SP100-T040S
		6.5 mm	10785SP100-T065S
		9 mm	10785SP100-T090S
Multi-Mode	SMA	3 mm	10785MP350-T030S
		4 mm	10785MP350-T040S
		6.5 mm	10785MP350-T065S
		9 mm	10785MP350-T090S
Multi-Mode	SMA	3 mm	10785MP450-T030S
		4 mm	10785MP450-T040S
		6.5 mm	10785MP450-T065S
		9 mm	10785MP450-T090S
	Single-Mode Multi-Mode	Single-Mode SMA Multi-Mode SMA	Single-Mode SMA 3 mm 4 mm Single-Mode SMA 6.5 mm 9 mm 9 mm Multi-Mode SMA 3 mm 4 mm Multi-Mode SMA 3 mm 6.5 mm Multi-Mode SMA 6.5 mm Multi-Mode SMA 6.5 mm

Part Numbering Schema

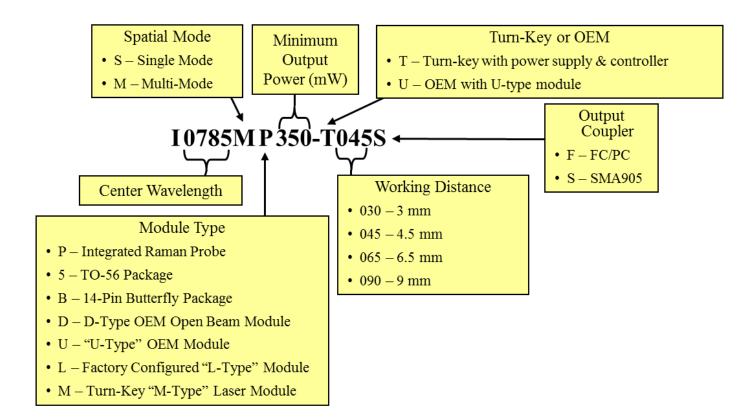





Innovative Photonic Solutions, Inc. 4250 U.S. Highway 1, Suite 1 Monmouth Junction, NJ 08852 Phone: (732) 355-9300 Fax (732) 355-9302 sales@innovativephotonics.com http://www.innovativephotonics.com

Mechanical Specifications Integrated Raman Probe

Probe Head

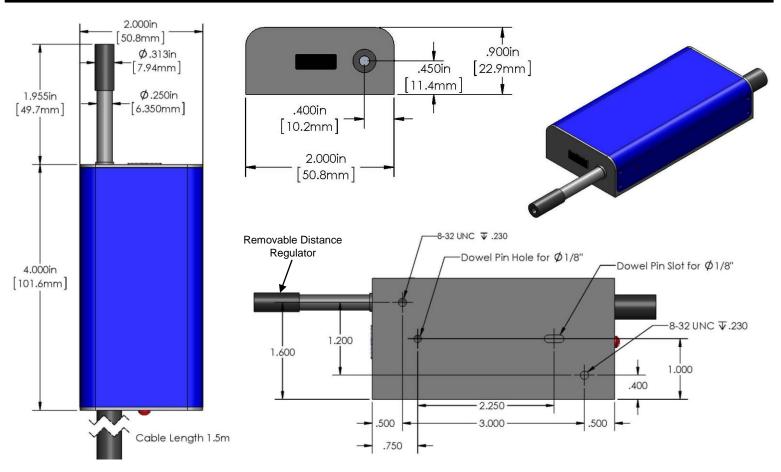

Control Module

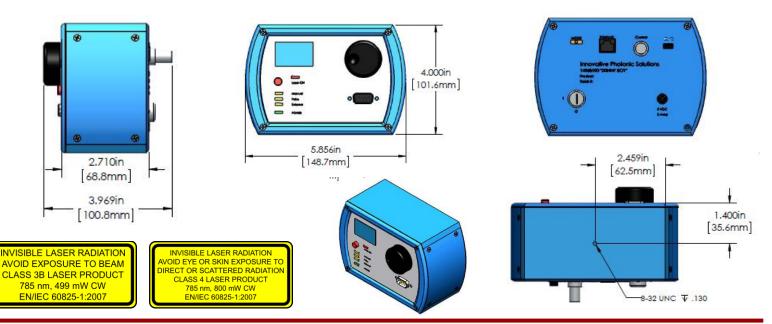

All data and statements contained herein are subject to change in accordance with Innovative Photonic Solution's policy of continual product improvement. No information contained herein is intended for use in connection with any contract except as may be first confirmed in writing by Innovative Photonic Solutions. The publication of information in this document does not imply freedom from patent or other rights of Innovative Photonic Solutions or others.